



Install Instructions for VC-100 Software V6.0

1. Click the Download link, to download the file to a temporary location on your computer.
2. Go to the temporary location on your computer, and unzip the file. (See bottom of this document for directions on how to unzip a file.)
3. Launch the setup.exe, if you haven't already done so, from your temporary location.
4. You will receive a few prompts from the program, and answer them to fit the install needs of your location.
5. When the installation has finished, a dialog box states that VC100WIN has completed successfully.
6. To run the program, go to Programs→VC100WIN→VC100 Version 6.0.
7. To send a shortcut to your desktop, go to Programs→VC100Win, and right-click on VC100 Version 6.0. Choose Send To→Desktop (create shortcut).

Instructions for .ZIP files

1. After downloading a .ZIP file, you must unzip it with [WinZip](#), [Pkunzip](#), or a similar program.
 - With WinZip, open the ".ZIP" file and select the "Install" function or icon. This extracts the files into a temporary directory and runs the "setup" program to start the installation. (The temporary files are then automatically deleted after the installation is done.)
 - With Pkunzip for Windows, open the ".ZIP" file and first extract the file into a temporary directory such as "c:\temp". Next, run the "setup" program to start the installation.
 - For a similar program, you must be familiar with the functionality of that program. We cannot cover all tools.
2. Run the "setup" program to install the Windows version.
3. Continue with the instructions above, at step number three.